

# Studying Structural Changes of Rod-like Nanoparticles Using Dynamic Light Scattering



Submitted by Viet Lam and Lynn Walker,

Department of Chemical Engineering, (Center for Complex Fluids Engineering), Carnegie Mellon University, Pittsburgh PA 15213

## Introduction

Polyelectrolyte-surfactant complexes are used in numerous applications ranging from detergents (for solubilization of oils) to more exotic applications like the molecular design of templates for nanomaterials. These self-assembled aggregates are extremely versatile, robust and often inexpensive. However, understanding the effect of the tenuous interactions that define the complexes (electrostatic-hydrophobic balance) is vital to the design and success of these materials in current and future applications.

A series of novel polyelectrolyte-surfactant complexes were studied in order to probe several candidates as potential pH sensitive materials. At neutral pH, the complexes are water soluble, rod-like nanorods. However, gelation and longer-range interactions are observed as the pH is dropped to low values (approximately pH 1). This is proposed to be a change in the complex structure, and dynamic light scattering (DLS) is used to determine changes in the overall complex size in dilute solutions. Currently, since the shape is not known, the translational diffusion coefficient that is measured in DLS is used to calculate an effective hydrodynamic diameter of the complex. The hydrodynamic diameter is measured as the pH of the solution is reduced from neutral to acidic conditions.

## Experimental

DLS measurements were made on a Malvern Zetasizer Nano S with a detection angle of  $173^\circ$  using a 4mW He-Ne laser operating at a wavelength of 633nm. The polyelectrolyte-surfactant complexes were synthesized at Carnegie Mellon University and freeze dried for storage. Samples are later reconstituted in filtered deionized water to prepare for measurement. The polyelectrolyte has a molecular weight of approximately 350K g/mol and the resulting polyelectrolyte-surfactant complex has a molecular weight on the order of 1,000,000 g/mol.

## Results and Discussion

Figure 1, shows the hydrodynamic diameter measurement results for a series of polyelectrolyte-surfactant complexes in solution at different values of pH. In this case, low concentrations were used in order to extract hydrodynamic data on individual particles without any concerns about interparticle interactions. A sharp increase in overall complex size is observed as the pH drops below 2, consistent with the observation of gelation at higher concentrations. This phenomenon is reversible and is not due to inter-complex interactions. This explanation was excluded by diluting the complexes at various values of pH


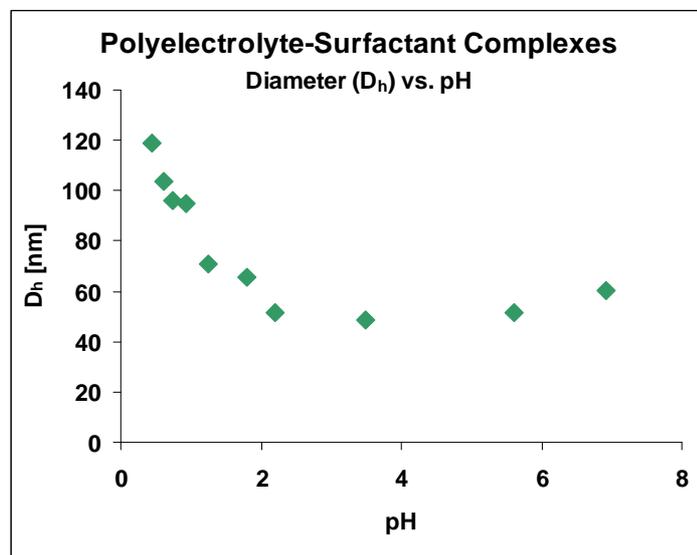


Figure 1: Hydrodynamic diameter measurements versus pH of polyelectrolyte-surfactant complexes in solution measured at 25°C

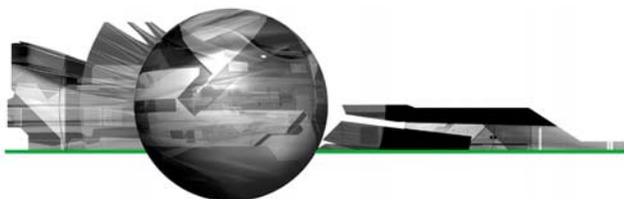


and finding no concentration dependence on size, which verifies that the change is actually a change in the complex size. This additional information suggests a change in structure which will be verified with other techniques (cryo-TEM or SANS).

The impact of temperature on the complex structure provides information about the relative strengths of electrostatic and hydrophobic interactions. Results shown in figure 2, indicate that the complex structure is not a function of temperature at neutral pH, but is a strong function of temperature at low pH. The structural changes in complex size as stated before are reversible and reproducible. These results are being used to develop a structural model for these materials.

The DLS results verify that the balance of the interactions controlling complex structure, change at low pH, indicated by the change in temperature dependence.

## References


1. Loxley, A. and B. Vincent, Equilibrium and kinetic aspects of the pH-dependent swelling of poly(2-vinylpyridine-co-styrene) microgels. *Colloid & Polymer Science*, 1997. 275(12): p. 1108-1114.
2. Yin, J., et al., pH-Induced Deswelling Kinetics of Sterically Stabilized Poly(2-vinylpyridine) Microgels Probed by Stopped-Flow Light Scattering. *Langmuir*, 2008.
3. Gerber, M.J., S.R. Kline, and L.M. Walker, Characterization of rodlike aggregates generated from a cationic surfactant and a polymerizable counterion. *Langmuir*, 2004. 20(20): p. 8510-8516.
4. Kuntz, D.M. and L.M. Walker, Solution behavior of rod-like polyelectrolyte-surfactant aggregates polymerized from wormlike micelles.

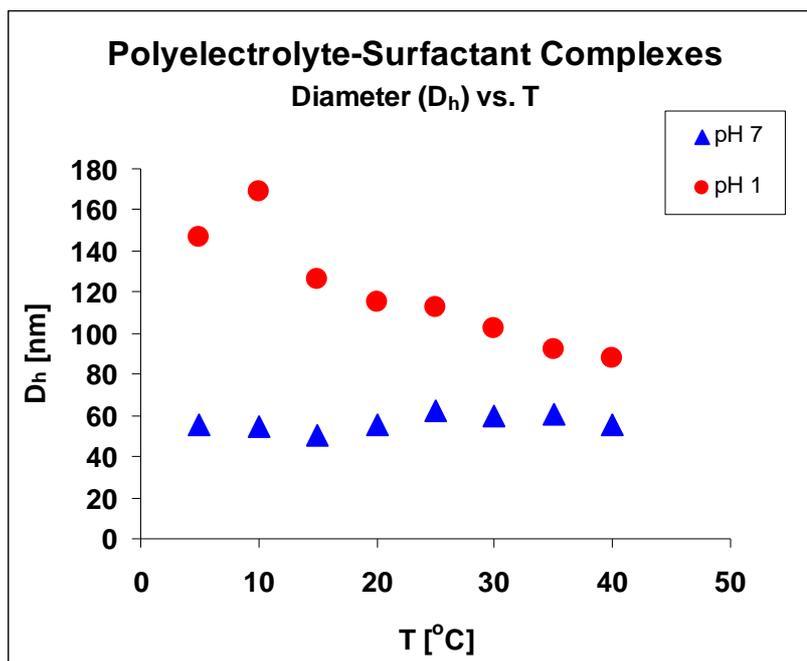

**Figure 2:** Hydrodynamic diameter measurements vs. temperature of polyelectrolyte-surfactant complexes under neutral (pH 7) and acidic (pH1) conditions

*Journal of Physical Chemistry B*, 2007. 111(23): p. 6417-6424.

5. Lam, V.D., Polyelectrolyte-surfactant complex – a pH induced transformation from rod-like to string-of-pearl structures, in preparation.

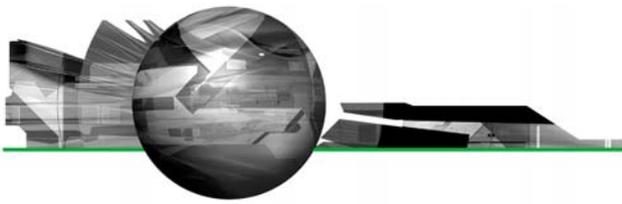

## Zetasizer Nano System

The Malvern Zetasizer Nano has temperature and time trend capabilities directly accessible in its software. In this particular study the Zetasizer Nano was used to measure the change in hydrodynamic diameter as a function of temperature. The Zetasizer Nano software is uniquely designed to make these measurements in an automated manner. By selecting a “trend measurement” it is easy to specify any temperature vs. size measurement profile from 2°C to 90°C.

Similarly, sample stability as a function of time can also be monitored. The same “trend” software will easily allow measurements of hydrodynamic diameter vs. time to be obtained. The addition of an automated multipurpose titrator, MPT-2, can also allow for automated pH or additives such as salts or surfactants or other processing aids to be studied as a function of size or zeta potential.

### Malvern Instruments Ltd

Enigma Business Park • Grovewood Road • Malvern • Worcestershire • UK • WR14 1XZ  
Tel: +44 (0)1684 892456 • Fax: +44 (0)1684 892789

### Malvern Instruments Worldwide

Sales and service centers in over 50 countries for details visit [www.malvern.com/contact](http://www.malvern.com/contact)

© Malvern Instruments Ltd 2009

[more information at www.malvern.com](http://www.malvern.com)